


LC-10ATvp

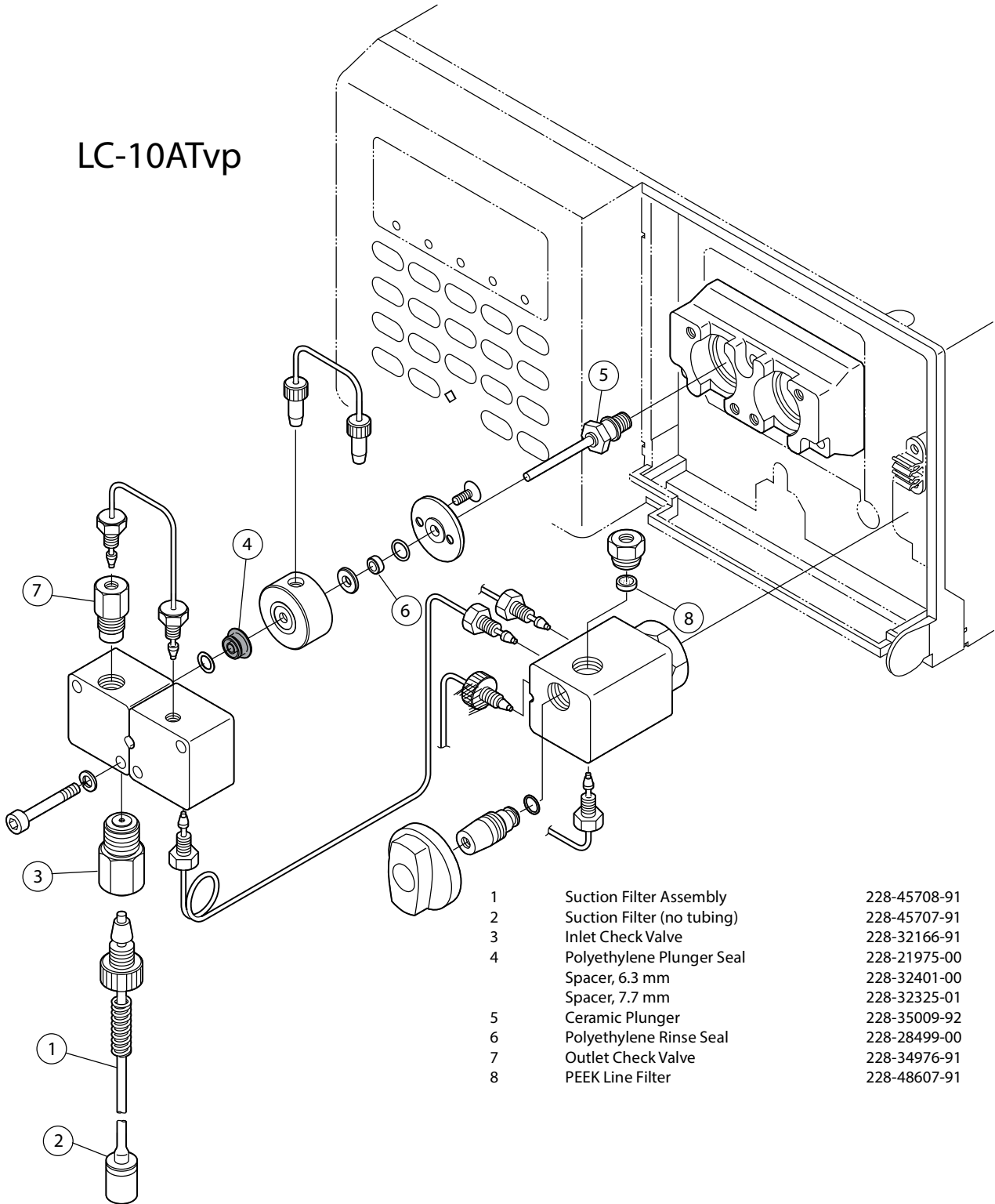

- | | | |
|---|----------------------------|--------------|
| 1 | Suction Filter Assembly | 228-45708-91 |
| 2 | Suction Filter (no tubing) | 228-45707-91 |
| 3 | Inlet Check Valve | 228-32166-91 |
| 4 | Polyethylene Plunger Seal | 228-21975-00 |
| | Spacer, 6.3 mm | 228-32401-00 |
| | Spacer, 7.7 mm | 228-32325-01 |
| 5 | Ceramic Plunger | 228-35009-92 |
| 6 | Polyethylene Rinse Seal | 228-28499-00 |
| 7 | Outlet Check Valve | 228-34976-91 |
| 8 | PEEK Line Filter | 228-48607-91 |