


# LC-20AD LC-20ADXR

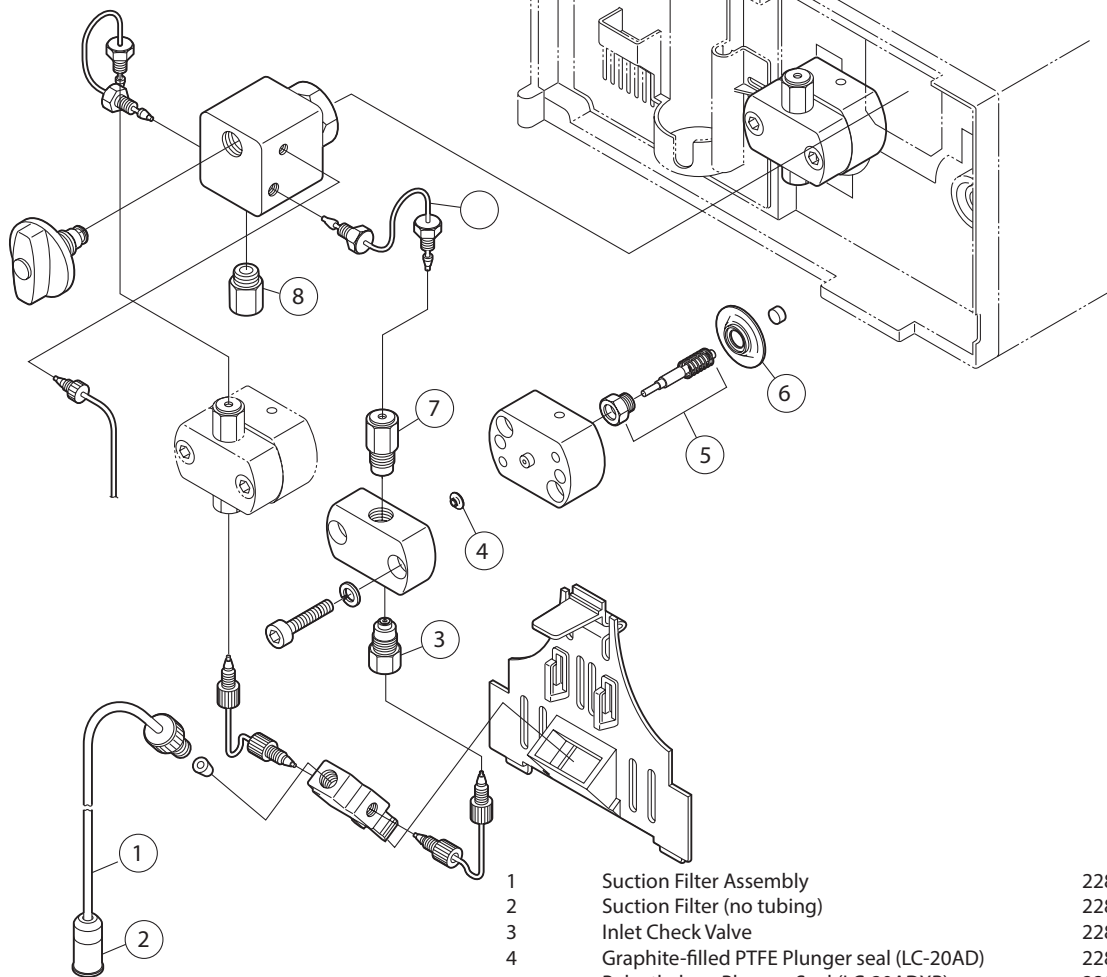

- |   |   |              |
|---|---|--------------|
| 1 | Suction Filter Assembly                     | 228-45708-91 |
| 2 | Suction Filter (no tubing)                  | 228-45707-91 |
| 3 | Inlet Check Valve                           | 228-48249-91 |
| 4 | Graphite-filled PTFE Plunger seal (LC-20AD) | 228-35146-00 |
|   | Polyethylene Plunger Seal (LC-20ADXR)       | 228-32628-91 |
| 5 | Sapphire Plunger                            | 228-35601-93 |
| 6 | Diaphragm (2 pack)                          | 228-32784-91 |
| 7 | Outlet Check Valve                          | 228-45705-91 |
| 8 | Line Filter                                 | 228-35871-96 |